(GB)

Biosart® 100 Monitor – Instructions for use Sartorius 100 ml Microbiological Monitor
Sterile, disposable 100 ml Microbiological Monitors with a 47 mm or 56 mm diameter membrane and large bottom port provides a faster flow rate and decreases testing time. Ideal for the detection of microorganisms in beverages, pharmaceuticals and water.

Materials included:

- 48 - Sterile, disposable microbiological monitors
- 48 - Bottom port plugs
- 1 - Large funnel adapter

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
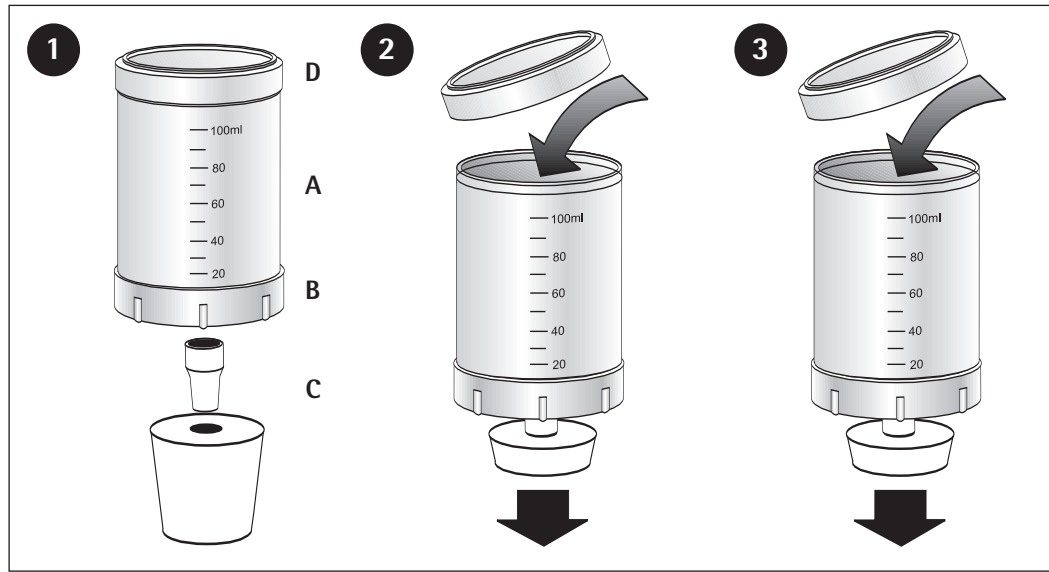
(GB)

1. Remove monitor from carton. Attach funnel adapter **C** to monitor base **B**. Securely insert the assembled unit to vacuum source.
2. Remove the lid **D**, add sample and replace the lid. Attention: if there is only a small volume to filter, pour sterile buffer solution (10 - 20 ml) into the monitor before pouring the sample. Apply vacuum and filter the solution.
3. Turn vacuum off and rinse the inside walls of the funnel with sterile buffer solution (20 - 30 ml). Apply vacuum. When filtration is complete, turn vacuum off. Release vacuum from the unit with either the vacuum release cap on a suction manifold or with a three-way stopcock when utilizing sidearm vacuum flask.
4. Remove the lid of the funnel and add your selected media ampoule to the membrane and pad from the top. Be careful not to touch the tip of the media ampoule to the membrane. Replace lid. Apply a one-second vacuum to draw the media down into the pad.
5. Holding funnel **A** firmly, twist off the monitor base **B**. Discard the funnel **A**, retain the funnel lid **D** and funnel adapter **C**.
6. Push the funnel lid **D** on to the monitor base **B**. Remove the funnel adapter **C** from the monitor and replace with a port cap **F**. Invert the monitor and incubate at appropriate temperature for selected media. Ensure that the horizontal position of the petri dish is exact so there is a homogenous distribution of the media.

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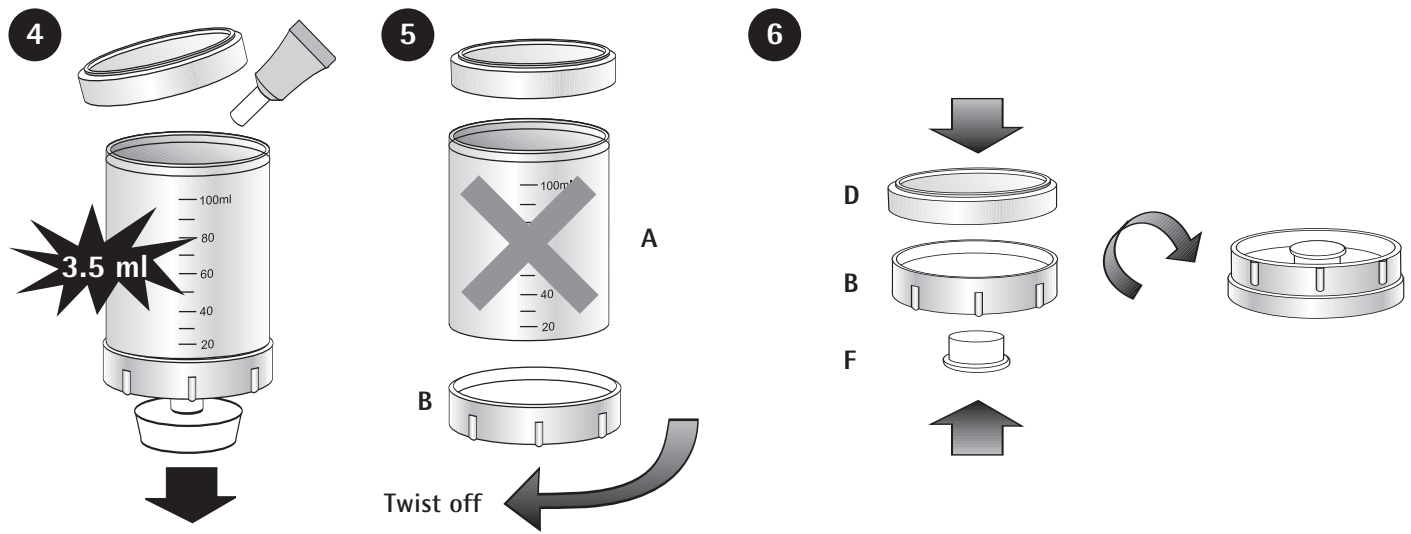Sales and Service Contacts

For further contacts, visit www.sartorius.com

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Phone +41.1.746.50.00
Fax +41.1.746.50.50

U.K.

Sartorius Ltd.
Longmead Business Park
Blenheim Road, Epsom
Surrey KT19 9QQ

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Phone +852.2774.2678
Fax +852.2766.3526

India

Sartorius India Private Ltd.
10, 6th Main, 3rd Phase Peenya
KIADB Industrial Area
Bangalore - 560 058

Phone +91.80.2839.1963/0461
Fax +91.80.2839.8262

Japan

Sartorius K.K.
KY Building., 8-11
Kita Shinagawa 1-chome
Shinagawa-ku
Tokyo 140-0001

Phone +81.3.3740.5407
Fax +81.3.3740.5406

Korea

Sartorius Korea
17-2 Jungja-Dong, Bundang-Gu
Sungnam, Gyunggi-Do
B-1023, Paragon
463-811, Korea

Phone +82.31.782.7011
Fax +82.31.782.7090

Malaysia

Sartorius (Malaysia) Sdn. Bhd.
Lot L3-E-3B, Enterprise 4
Technology Park Malaysia
Bukit Jalil
57000 Kuala Lumpur

Phone +60.3.8996.0622
Fax +60.3.8996.0755

Sartorius Singapore Pte. Ltd.
10, Science Park Road, The Alpha
#02-25, Singapore Science Park 2
Singapore 117684

Phone +65.6872.3966
Fax +65.6778.2494

Australia

Sartorius Australia Pty. Ltd.
Unit 17/104 Ferntree Gully Road
Waverley Business Park
East Oakleigh, Victoria 3166

Phone +61.3.9590.8800
Fax +61.3.9590.8828

Biosart® 100 Monitor

Instruction Manual