



■ Introducing the Labplas SANI-STICK



- Ideal for surface sampling
- Allows to reach hard-to-reach areas
- Easier handling
- One time use
- Made to avoid sponge wedging
- Plastic sponge handle with a strong grip
- Comes with a 2.5 mil.in sterile sampling bag; 4.5 x 9 in
- **TWIRL'EM** bag with leak proof closure system
- Sterilized by gamma irradiation
- 100 units per case (10 packs of 10 units)
- **Choice of:**
 - Neutralizing agent
 - Neutralizing agent DE



Sani-stick user guide

STEP 1	STEP 2	STEP 3	STEP 4	STEP 5
				
Remove the tear-off strip and carefully take the sani-stick by the handle.	The wet sponge is already on the handle; swipe it on the determined surface.	Press on the bottom of the handle to gently release the sponge in the sterile sampling bag.	Twirl the bag three times and close it with the tabs.	Identify the sample on the write-on area with a permanent marker.

Catalog No.	Description	mil.in	micr.	in x in	mm x mm	oz	ml	Packaging
KSS61310NE	Sponge on handle with neutralizing buffer	2.5	63	4.5 x 9	114 x 229	15	450	Cs/100 (10 x 10)
KSS61310DE	Sponge on handle with DE neutralizing buffer	2.5	63	4.5 x 9	114 x 229	15	450	Cs/100 (10 x 10)

■ STERILITY CERTIFICATE IN EVERY BOX/FDA GRADE PLASTIC

CORPORATE HEADQUARTERS

1430 West McCoy Lane
Santa Maria, California
United States 93455

CUSTOMER SERVICE

Phone:
805-346-2766 or 800-266-2222
Fax:
805-346-2760

